


pi-pump User Instructions

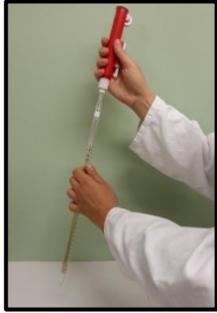

1. Carefully insert your pipette into the cone while applying light pressure.




2. Put top of the pipette into the liquid.



3. Turn the knurled wheel down to suck up the liquid.



4. Pressing the lever will release all the liquid instantly.

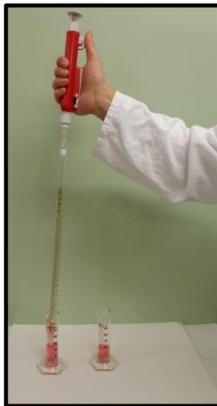

5. To release the liquid more precisely, i.e. line by line, turn back the wheel.



6. Pressing down the piston can release the liquid as well.



7. Once you've finished your work, pull the pipette gently to remove it from the cone.

Be careful when inserting and removing the pipette.
Extreme care should be taken when inserting disposable glass pipettes, as they can break.

Glasfirn Giessen

Inh. Sabine Körber e.K.
Kirchstraße 34
D-35396 Gießen-Wieseck
Tel.: +49 (0)641-5832
Fax: +49 (0)641-5833
E-Mail: info@glasfirn.de
www.glasfirn.de

The pi-pump is available in four sizes (colour-coded):
For pipettes of up to:
- 25ml, red, order no. 2524
- 10ml, green, order no. 2520
- 2ml, blue, order no. 2516
- 0.2ml, yellow, order no. 2512 (without quick release)